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The apparatus is used for the verification and quantitative determination of the linear expansion occurring due to temperature changes on tubes of different material and different lengths.

Bibliography:

Central Card File Physics Demonstration,  
Heat (598 631, in German)

#### 1 Safety notes

Hot water vapour!  
Hot tubes! Use protective cloth when changing tubes!

#### 2 Scope of supply, technical data

- ① Basic apparatus, 73 cm x 5 cm x 3 cm
- ② Screw for stand rod (for mounting the expansion apparatus according to Figs. 3 and 4)
- ③ Screw for fitting the tube into the fixed bearing ④
- ④ Fixed bearing for the fixed tube end, tube attached by means of the screw ⑤ at 600, 400 or 200 mark of the basic apparatus so as to obtain effective tube lengths of 600 mm, 400 mm or 200 mm
- ⑤ Screw to mount bearing ④ onto the basic apparatus
- ⑥ Projectable scale, 14 cm x 20 cm, 0 to 1 cm, 0.05-mm graduation
- ⑦ Pointer connected by rack and pinion with guide bearing ⑧
- ⑧ Guide bearing for movable tube end
- ⑨ Spring clips to secure the tubing
- ⑩ Tubes of steel, brass and glass, each 65 cm long, 7 mm O. D., with annular grooves mating the knurled-head screw ③ of the fixed bearing ④ and with hose nozzles to connect tubing of 7 mm I. D.

(on delivery of the apparatus, two tubes are fixed below the basic apparatus ① by means of clamps)

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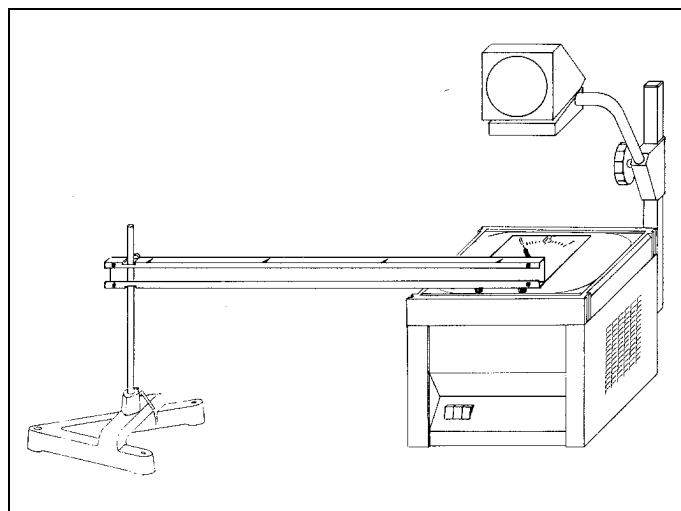
Fig. 3

### 3 Operation

#### 3.1 Additionally required equipment for heating the tubes:

*Heating by water steam (see Fig. 2)*

1 Tripod	666 683
1 Wire gauze, heat resistant	666 685
1 Bunsen burner, e.g. Bunsen burner for butane gas	666 711/712
1 Erlenmeyer flask	664 233
1 Rubber stopper with 1 hole	667 265
1 Glass tube, approx. 8 mm O. D. e.g.	665 204
Rubber tubing	307 65

or

1 Steam generator	303 28
-------------------	--------

or

1 Lid with heater	590 50
1 Aluminium beaker	590 05
1 Low-voltage power supply, e.g.	591 09
1 Beaker as collector, e.g. from	664 103

#### 3.2 Modes of reading and corresponding set-up:

*Direct reading of the changes in length on the scale of the expansion apparatus (Fig. 2)*

Recommended equipment:

Laboratory stand	300 75
------------------	--------

Additional equipment required for projecting the scale:

*For shadow projection:*

Lamp housing with lamp, 6 V, 30 W	450 60/51
Transformer, 6 V, 30 W, e.g.	562 73
Saddle base	300 11

*Reading from the dial gauge (Fig. 4):*

(especially recommended where high measuring accuracy is required)

Dial gauge	361 15
Holder for dial gauge	361 36
Laboratory stand	300 75
Stand material	

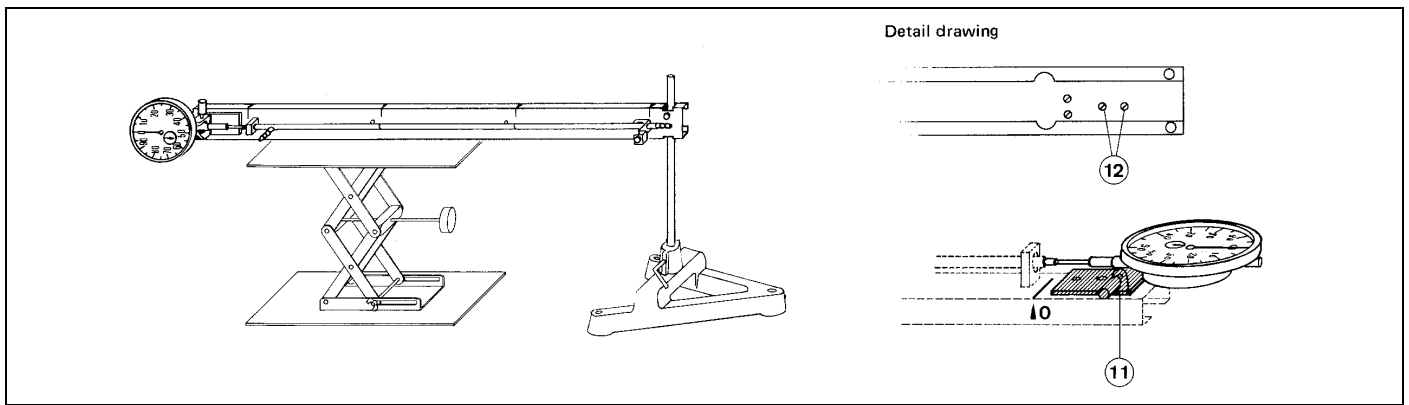
Fig. 4

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To exchange the projectable scale ⑥ with the dial gauge, proceed as follows (Fig. 4):

Remove scale and place the instrument on its side (see detail drawing). Demount the contact pin with rack and pinion gear and pointer by loosening the screws ⑫, and replace by dial gauge ⑪. Clamp the dial gauge in the holder ⑪.

### 3.3 Preparing and conducting the experiment

Fasten the fixed bearing ④ at mark 600, 400 or 200.

Slide the open tube end into the fixed bearing and the closed tube end into the guide bearing ⑧. Fix the tube by tightening screw ③.

*Important:* The screw should engage into the annular groove of the tube.

After connecting the rubber tubings, set the pointer ⑦ to zero according to Fig. 2.

Place a collecting vessel below the tube as shown in Fig. 2.

## 4 Storage

It is advisable to clamp the glass tube below the basic apparatus in order to avoid damage (e.g. by pressure or shock).