
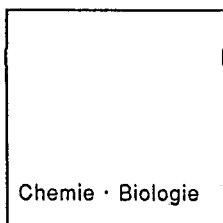
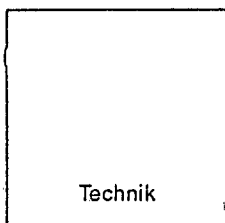
Physik

Chemie · Biologie

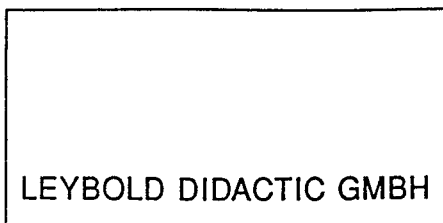Technik

LEYBOLD DIDACTIC GMBH

12/91-kem-

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The adhesive magnetic board serves as base plate for demonstration experiments with magnetically adhesive units, .e.g.

- in geometrical optics with unit 463 76 or the collection 463 70/71,
- in mechanics with the magnetic bases 301 330 or the collection 343 001,

and as a background plate for experiment set-ups in the demonstration panel frame (301 300).

#### 1 Safety instructions

- Use only water-soluble markers for labelling.
- Clean the board with water only or, if necessary, water with dishwashing liquid added. If necessary, clean residue with alcohol (spirits). Do not use detergents.
- In order to avoid scratching the board surface, take care that the magnetic surfaces of the magnetically adhesive units are free of dirt, particularly metallic particles.

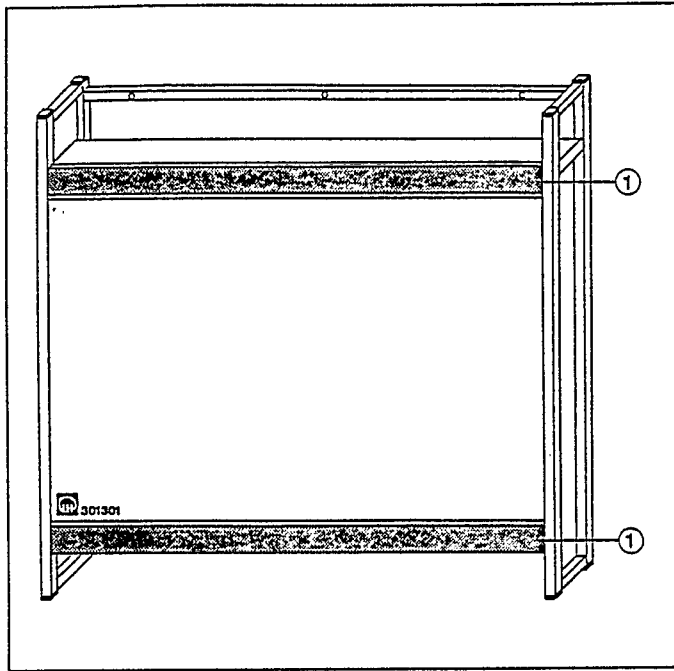
#### 2 Technical data

Ferromagnetic plate, with painted imprint

Working surface: 92.5 cm x 59 cm

Imprints: rastered grid, crosspoints separated by a distance of 5 cm  
Angular scale, 360°/10°, 20 cm Ø

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Fig.1 Magnetic board in experiment frame  
① Profile rails

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#### 3.1 Experiment frame (i.e. 301 300), cf. fig. 1

Mount the magnetic board in the profile rails (cf. fig. 1). To do this, first fit the top edge of the board into the upper profile rail, then fit the bottom edge into the lower profile rail.

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Fig.2 Mounting in assembly frame  
① Slotted block

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**3.2 Demonstration assembly frame (598 07) with pair of tracks (598 09/10), cf. fig. 2**

The slotted blocks (available separately, 200 69 833, includes 4 units) enable the adaption of tracks (598 09 or 598 10) to the demonstration assembly frame (598 07). The slotted blocks are screwed on to the magnetic board.

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Fig.3 Assembly with stand equipment  
① Metal plate

**3.3 Stand equipment, cf. fig. 3**

4 Leybold multiclamps (301 01) recommended, each with metal plate (200 65 559) can be screwed on for setting up to stand equipment.