05/03-W97-Hund

Installationsanleitung

Pocket-CASSY® (524 006)

1 Installation

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Pocket-CASSY® (524 006)

1 Installation

Pocket-CASSY is connected to a PC via a USB port using the enclosed USB cable. For this a 500 mA port is required. Usually all built-in ports as well as ports in USB hubs with external voltage supply (e.g. monitors, but not keyboards) are of this type.

When Pocket-CASSY is connected for the first time, it is identified by the operating system automatically and the necessary drivers are installed. Under Windows 98 the driver installation has still to be done manually. To do this, confirm the dialog windows that appear on the screen with "Continue" and finally insert the Windows installation CD ROM. After the drivers have been installed, Pocket-CASSY is found automatically by the CASSY Lab software (524 200).

CASSY Lab is installed automatically by inserting the CASSY Lab CD ROM or manually by starting the file SETUP.EXE and following the instructions on the screen. The software is configured under „Start → Programs → CASSY Lab“.

2 Technical data

Please refer to the CASSY Lab online documentation for the technical data.

In conjunction with Pocket-CASSY, CASSY Lab runs on Windows 98/ Me/ 2000/ XP systems with Pentium III, 300 MHz or higher. Several Pocket-CASSYs can be operated simultaneously on different USB ports. However, CASSY Lab cannot be used in conjunction with Pocket-CASSY and other CASSY modules (e.g. Sensor-CASSY) at the same time.

3 Online documentation

Please refer to the CASSY Lab online documentation for the program description of CASSY Lab and for the description of the different CASSY modules. In this documentation, numerous experiment examples are described as well.

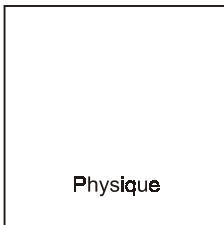
Safety notes

At the present state of software development it is not possible to guarantee that a product is faultless. Therefore Leybold Didactic GmbH does not take responsibility for any damage resulting from the installation or use of Pocket-CASSY and CASSY Lab.


Electrical safety:

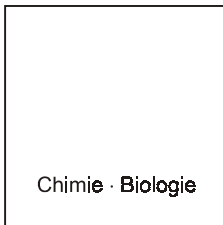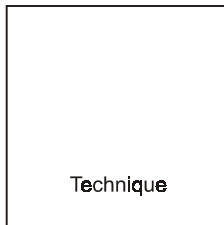
- Do not apply voltages over 30 V to sensors and sensor boxes which are connected to Pocket-CASSY.
- Pay attention to possible ground connections between sensors and the computer.

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